

Manual

Proteinase K

Enzyme for digesting proteins in biological samples. Concentration 20 mg/ml. Activity ≥ 30 U/mg.

catalog #	size
1019-20	1 ml
1019-20-5	5 x 1 ml

For research use only.

Guarantee

A&A Biotechnology provides guarantee on this product.

The company does not guarantee correct performance of this kit in the event of:

- not adhering to the supplied protocol
- use of not recommended equipment or materials
- use of other reagents than recommended or which are not a component of the product
- use of expired or improperly stored product or its components

Description

Proteinase K (EC 3.4.21.64) is a broad-spectrum serine protease. The predominant site of cleavage is the peptide bond adjacent to the carboxyl group of aliphatic and aromatic amino acids with blocked alpha amino groups.

Proteinase K is inactivated by diisopropyl fluorophosphate (DFP) or phenyl methane sulfonyl fluoride (PMSF).

It's a recombinant enzyme cloned from fungus *Engyodontium album* and produced in *Pichia pastoris*.

Application

- extremely effective proteolytic degradation of biological material
- inactivation of endogenous DNAses and RNAses during nucleic acid purification

Contents

	1019-20	1019-20-5	storage
proteinase K, recombinant	1 ml	5 x 1 ml	4-8 °C
storage buffer: 20 mM Tris, pH 7.5, 1 mM CaCl ₂ , 0.02% sodium azide, 50% glicerol (v/v)			

Unit definition

1 U of proteinase K releases 1 µmole of Folin positive amino acid in 10 min at 37 °C, pH 7.5, using denatured hemoglobin as a substrate.

References

1. Burkiewicz A., Dąbrowski S., Barski P., *Polska rekombinowana Proteinaza K*, (2007) *Postępy Biochemii* 53(4): 327-328
2. Ebeling W., Hennrich N., Klockow M., Metz H., Orth H.D., Lang H., (1974) *Eur. J. Biochem.* 47(1):91-97
3. Hilz H., Wiegers U., Adamietz P., (1975) *Eur. J. Biochem.* 56(1):03-08
4. Betzel C., Sigh T.P., Visanji M., et al., (July 1993) *J. Biol. Chem.* 268(21):15854-15858

Safety information



DANGER

Proteinase K

H315 Causes skin irritation.

H319 Causes serious eye irritation.

H334 May cause allergy or asthma symptoms or breathing difficulties if inhaled.

H335 May cause respiratory irritation.

P261 Avoid breathing dust.

P305+P351+P338 If in eyes: rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

P342+P311 If experiencing respiratory symptoms call a Poison Center or doctor/physician.



A&A BIOTECHNOLOGY

innovating life science

A&A Biotechnology, ul. Strzelca 40, 80-299 Gdańsk, Poland
phone: +48 883 323 761, +48 600 776 268
info@aabiotech.com, www.aabiotech.com

version 2024-1

