

Manual

MagnifiQ™ 1 Pathogen instant kit

Kit for automated, magnetic isolation of DNA and RNA of pathogens in the strip format. Contains ready-to-use, reagent-filled stripes and all necessary consumables.

catalog #	size	compatible devices *
607A-1V-32	32 isolations	Auto-Pure Mini
607A-1V-160	160 isolations	Auto-Pure Mini

* Compatible devices

The kit has been tested with specific Allsheng brand isolation devices. This does not preclude it from working with other devices. If your device is not listed, please contact us at info@aabiotech.com.

For research use only.

Guarantee

A&A Biotechnology provides a guarantee on this product.

The company does not guarantee the correct performance of this kit in the event of:

- not adhering to the supplied protocol
- use of not recommended equipment or materials
- use of other reagents than recommended or which are not a component of the product
- use of expired or improperly stored product or its components.

Table of Contents

Advantages	3
Specification	3
Description	3
Contents	3
Additional equipment and reagents	4
Necessary	4
Optional	4
Material preparation	4
Swabs with transport medium	4
Dry swabs	4
Protokol	5
Safety information	7

Advantages

- Requires only a few minutes of manual work while adding the samples. The rest of the procedure is carried out in an automatic extraction device.
- It does not require the initial preparation of buffers. Just add samples to the strip and after about 30 min you get extracted material.
- Universal extraction of both DNA and RNA.

Specification

protocol time	~30 min
sample type	swabs
sample size	up to 400 µl
elution volume	100 µl
elution solution	Tris buffer
binding capacity	30 µg DNA and RNA
downstream applications	sequencing, qPCR, RT-PCR

Description

MagnifiQ™ 1 Pathogen instant kit is designed for RNA and DNA isolation from viruses and Gram(-) bacteria. The isolated material is perfect for further analyzes and tests by qPCR and RT-PCR methods and for sequencing.

The MagnifiQ™ product series is based on the automated isolation of nucleic acids with use of magnetic beads. This method significantly shortens working time and reduces risk of mistake in comparison to manual methods.

Contents

component	607A-1V-32		607A-1V-160		storage
	quantity	cat #	quantity	cat #	
XS-G - extraction strip	4 x 8 pcs	K-S1V22XG	20 x 8 pcs	K-S1V22XG	15–25 °C
Proteinase K	1,1 ml	K-PRK-11A	4 x 1,1 ml	K-PRK-11A	4–8 °C*
tip comb 8	8 x 2 szt.	K-C8U-2	40 x 2 szt.	K-C8U-2	15–25 °C

* Proteinase K can be stored at 15–25 °C for up to 12 months.

Additional equipment and reagents

Necessary

- pipette
- pipette tips
- 2 ml Eppendorf tubes (sample lysis)

Optional

- vortex
- sterile water, Tris buffer, PBS buffer

Material preparation

Swabs with transport medium

No additional material preparation is required.

Dry swabs

1. Break or cut off part of the swab with the collected sample and place it in a 2 ml Eppendorf tube (not included).

Note. The portion of the swab with the collected sample should fit completely into the tube.

2. Add 500 µl of sterile water, Tris buffer or PBS buffer.


Note. Part of the swab with the sample should be completely immersed in the buffer.

3. Vortex for 20 s..

4. Follow point 1. [of the Protocol](#).

Protokol

1. Place **XS-G** stripes in the rack.

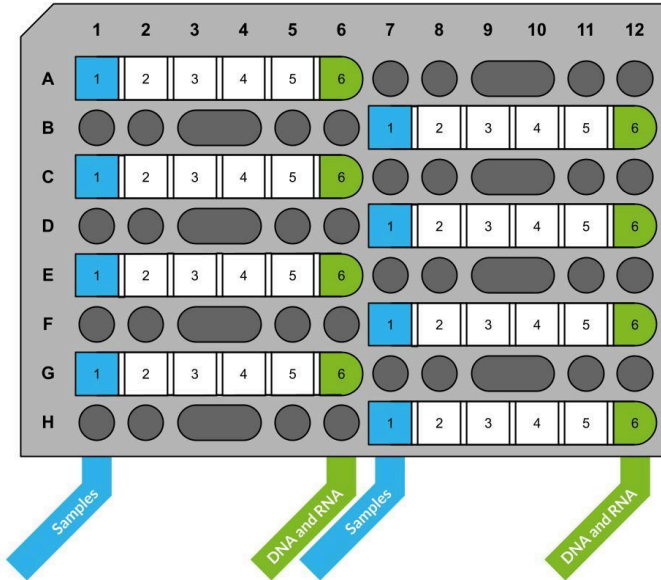

2. Remove the foil from the **XS-G** stripes.
3. Add **200-400 µl** of samples to the well **1** (first from the left) on the **XS-G** strip.
Note. The wells are numbered on the side of the strip.
Optional. Add 20 µl of Proteinase K to the wells 1 on the XS-G strip.
4. Place the rack in the extraction device.
5. Place the appropriate number of **tip combs 8** in the extraction device.

6. Run the program according to the table below:

Step	Well	Name	Mix Time (min)	Magnet (s)	Wait Time (min)	Volume (µl)	Mix Speed (1-10)	Temp. (°C)
1	1	LYSIS	5,0	0	0,0	900	4	50
2	5	BEADS	0,5	30	0,0	600	5	OFF
3	1	BIND	5,0	60	0,0	900	6	OFF
4	2	WASH1	2,0	30	0,0	600	6	OFF
5	3	WASH2	1,0	30	0,0	600	6	OFF
6	4	WASH3	1,0	30	5,0	600	6	OFF
7	6	ELUTION	6,0	60	0,0	100	8	65
8	3	DROP	1,0	0	0,0	600	5	OFF

7. After the program is over, remove the rack from the extraction device and transfer the purified DNA and RNA located in well **6** (first from the right) on the **XS-G** strip into sterile tubes (not included).

Note. Store extracted material at 4 °C for DNA or -70 °C for RNA.

Safety information



DANGER

Proteinase K

H315 Causes skin irritation.

H319 Causes serious eye irritation.

H334 May cause allergy or asthma symptoms or breathing difficulties if inhaled.


H335 May cause respiratory irritation.

P261 Avoid breathing dust.

P305+P351+P338 If in eyes: rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

P342+P311 If experiencing respiratory symptoms call a Poison Center or doctor/physician.

XS-G extraction strip

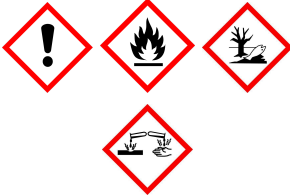

DANGER

H225 Highly flammable liquid and vapor.

H302+H312+H332 Harmful if swallowed, in contact with skin or if inhaled.

H314 Causes severe skin burns and eye damage.

H412 Harmful to aquatic life with long lasting effects.

P210 Keep away from heat, hot surfaces, sparks, open flames and other ignition sources. No smoking.

P273 Avoid release to the environment.

P280 Wear protective gloves/protective clothing/eye protection/face protection/hearing protection.

P301+P312+P330 If swallowed: Call a poison center/doctor/ if you feel unwell.

P303+P361+P353 If on skin (or hair): Take off immediately all contaminated clothing. Rinse skin with water or shower.

P304+P340 If inhaled: Remove person to fresh air and keep comfortable for breathing.

P305+P351+P338 If in eyes: rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.



A&A BIOTECHNOLOGY

innovating life science

A&A Biotechnology, Strzelca 40, 80-299 Gdańsk
tel. 883 323 761, 600 776 268
info@aabiotech.com, www.aabiotech.com

version 2024-1

